
Manual for HM-100S Overhead Stirrer



Thanks so much for your order of our overhead stirrer. In order to use freely the stirrer, please read this manual before using. Thanks.

I. Product illustration

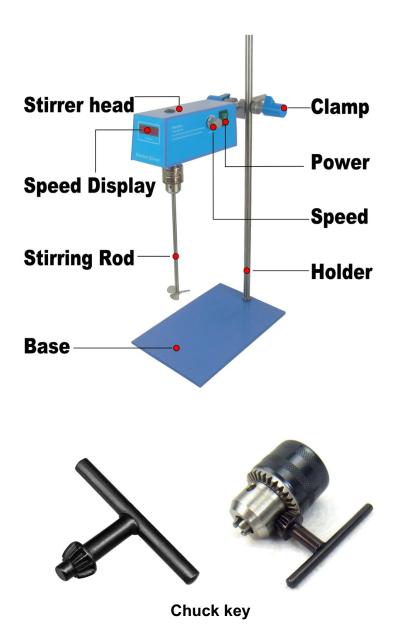

II. Installation

- 1. Use the holder to the base to fix tightly.
- 2. Installed Clamp to the holder and keep the one port is up so that we can fix the stirrer head later.
- 3. Take the stirrer head from the package and use the chuck Key to fix the stirring rod tightly.
- 4. Please fix the stirrer head to the clamp.
- 5. Please connect to power.
- 6. Turn on the Power and use the speed button to adjust the speed you need, the LCD screen will be showed.

III. After sales service

During use, any questions, please contact us. We will help you to solve it ASAP.

Contact information:

Shenzhen Haocheng Instrument Co.,Ltd

Attn: Steven

Email: hcinstrument@163.com

Mob./What's App: +86-134-2513-4344